



M3:10AC™ Supplemented Culture Media (with 10% Fetal Bovine Serum and Antibiotics)

Product Name: M3:10AC™
Product Codes: M310AC-100, M310AC-500
Product Use: Serum-free, highly enriched liquid culture medium with various growth factors and supplements; for human stem cells and other progenitor and fastidious cell types; member of M3™ media family
Features: GMP quality; sterile; USP grade materials; contains antibiotics

Manufacturer: INCELL Corporation LLC
Address: 12734 Cimarron Path
City/State/Zip: San Antonio, TX 78249
Phone: 800.364.1765; 210.877.0100
Email: info@incell.com
FAX: 210.877.0200

Distributor: Charter Medical
Address: 3948-A Westpoint Boulevard
City/State/Zip: Winston-Salem, NC 27103
Phone: 866.458.3116 (Toll Free)
Email: info@chartermedical.com
FAX: 336.714.4241

General Description

M3:10AC is a proprietary, multi-factor, supplemented high performance enriched culture media in the M3 Media Family (Table 1). M3:10AC™ contains 10% fetal bovine serum (FBS) and antibiotics, but no phenol red. It supports the in vitro growth of human stem and progenitor cells, many fastidious cell types, including primary human tumor cells and tissues. Its counterpart formula with phenol red is M3:10A™.

Formulation and Packaging

GMP manufactured M3:10AC™ is an enriched, supplemented basal media that contains salts, amino acids, peptides, multiple growth-promoting supplements, and antibiotics (clindamycin, gentamicin, amphotericin B). M3:10AC™ media are packaged in 100 mL and 500 mL bottles. Additional serum or other factors can be added per the needs of the tissue or cell type, and the intended use. The M3™ media family members are shown in Table 1. M3:BaseAC™ is the enriched formula to which the 10% v/v FBS is added to make the M3:10AC™ media. M3:DC™ (no phenol red) or M3:D™ are commonly used as "control", "holding" or "shift-down" media in studies where M3™ media are being used for growth or in vitro testing, and a control group requires absence of growth stimulating factors.

Use and Methods

As summarized in Tables 2 and 3, respectively focused on humans and animals, an extensive experience with many types of tissues and cells released for growth in culture have successfully used M3:10™ media. A variety of methods to support cell monolayers and/or suspension cell growth have been used with suitable M3™ supplemented media. Cell substrates have included standard cell culture plastics, complex biomatrices, meshes and engineered scaffolds.

Specifications

Visual
pH (USP <791>)
Osmolality (USP<785>)
Sterility: SC (USP <71>)
Sterility: FTG (USP <71>)
Mycoplasma (USP <63>)
Endotoxin (USP <85>)
Expiration

Acceptance Criteria

Clear, colorless
6.9 to 7.7
340 mmol/kg to 400 mmol/kg
No microbial growth
No microbial growth
None detected
<2.5 EU/mL
18 months from date of manufacture

Table 1. M3™ Media Family: Media Designations and Descriptions

Table with 4 columns: Media Designation, Item Number, Antibiotics, Phenol Red. Rows include M3:BaseA™, M3:BaseAC™, M3:BaseF™, M3:BaseFC™, M3:10A™, M3:10AC™, M3:20AC™, M3:D™, and M3:DC™.

The media suffix designations of "F", "A" and "C" refer, respectively, to "Free (F) of antibiotics", contains "Antibiotics" (A: gentamicin (50 µg/mL); clindamycin (6.5 µg/mL) and/ or Amphotericin B (2.5 µg/mL); are in the media, and Clear (C; no phenol red). Numbers after media designations are percent (v/v) serum in formula. Sera can be added to media at preferred concentrations.



M3:10AC™

Complete Culture Medium with 10% FBS and Antibiotics



## Manufacturing

M3:10AC™ is manufactured by sterile 0.22 µm filtration and packaging, using cGMP standards in an ISO Class 7 clean room and ISO Class 5 biosafety cabinet and using USP Standards for QC testing. Raw materials are pre-tested and the final product is checked for endotoxin, sterility (bacteria, fungi, and mycoplasma) and other quality specifications and criteria prior to release and preparation of the Certificate of Analysis.

**Table 2. Human Tissues and Cells in M3™ Media for Oncology and Regenerative Medicine Clinical and Research Applications**

| Human Tissues                             | Primary Culture Cell Types or Cell Lines [Media]  |
|---|---|
| Adipose (Fat)                             | Mesenchymal Stem Cells; Stromal vascular fraction regenerative cells [M3:10™]; adipose cells [M3:30™]   |
| Bone Marrow; Bone, Cartilage, Adipocytes  | Hematopoietic and mesenchymal stem cells; various types of renewable progenitor cells; Endothelial cells; entire population [M3:20™]; subsets of cells in other M3™ media; induced bone and cartilage and adipocyte outgrowth and/or induction of differentiation |
| Colon                                     | Primary epithelial and/or mesenchymal support cells [M3:2™] [M3:10™] and INCELL Cell Line NCM460 [M3:10™]   |
| Gastrointestinal                          | Primary epithelial cells and/or mesenchymal support cells; [M3:2™] [M3:5™] [M3:10™]   |
| Kidney                                    | Primary epithelial cells and/or mesenchymal support cells; [M3:2™] [M3:5™] [M3:10™]   |
| Liver                                     | Primary epithelial cells and/or mesenchymal support cells; [M3:2™] [M3:5™] [M3:10™]   |
| Muscles (Heart; Peripheral; Smooth)       | Pericytes; Mesenchymal or Stromal Stem Cells; regenerative cells [M3:10™]   |
| Nucleus pulposus (NP) Intervertebral Disc | NP stem cells; annulus chondrocytes and mesenchymal stem cells; various types progenitor cells [M3:10™]; etc.   |
| Pancreas                                  | Pancreatic islet beta and acinar cells [M3:5™]; etc.  |
| Peripheral or apheresis blood             | Circulating or mesenchymal cells; endothelial cells [M3:20™]; subsets of cells in other M3™ media formulas  |
| Placenta                                  | Hematopoietic, endothelial and mesenchymal stem cells; trophoblasts; syncytiotrophoblasts; various renewable progenitor cells; [M3:10™]; others   |
| Skin (adult; foreskin)                    | Epidermal keratinocytes co-cultures; Dermal Fibroblasts; Mesenchymal cells [M3:10™]; others   |
| Tumors; various                           | Epithelial, mesenchymal, lymphoid; [M3:10™]; others   |
| Umbilical Cord                            | Hematopoietic, endothelial and mesenchymal cells; various types of renewable progenitor cells; [M3:10™]; etc.   |

**Table 3. M3™ Complete, Supplemented Media Have Been Used for Animal Tissues and Cells Cultured for Biomedical Research**

| Cells and Tissues Derived from Adult, Newborn and/or Fetal Sources |  |
|--|--|
| Animal Tissues   | Species & Cultured Cell Types or Cell Lines [Media]  |
| Adipose (Fat)  | Rat, mouse, hamster, rabbit; Mesenchymal Stem Cells; Stromal vascular fraction regenerative cells [M3:10™]; adipose cells [M3:30™]*  |
| Bone Marrow  | Rat, mouse, hamster, rabbit; Hematopoietic and mesenchymal stem cells; various types of renewable progenitor cells; Endothelial cells; entire population [M3:20™]; subsets of cells in other M3™ media |
| Brain & Neural (Spinal)  | Rat, mouse: Progenitors; induced differentiation; [M3:5™] [M3:10™]   |
| Colon; Gastrointestinal  | Rat, mouse, hamster; Primary epithelial and/or mesenchymal support cells [M3:2™]; [M3:5™]; [M3:10™] or complex tissues in organ-like cultures  |
| Kidney   | Rat, mouse, hamster; Primary epithelial cells and/or mesenchymal support cells; [M3:2™]; [M3:5™]; [M3:10™]   |
| Liver  | Rat, mouse, hamster; Primary epithelial cells and/or mesenchymal support cells; [M3:2™]; [M3:5™]; [M3:10™]   |
| Muscles (Peripheral; Heart; Smooth)                                | Rat, mouse, hamster, rabbit; Pericytes; Mesenchymal or Stromal Stem Cells; regenerative cells [M3:10™]   |
| Pancreas; Other Neuroendocrine                                     | Rat, mouse, hamster; pancreatic islet beta and acinar cells [M3:5™]; other organs (e.g., adrenal)  |
| Peripheral or apheresis blood                                      | Rat, mouse, hamster, rabbit; Circulating or mesenchymal cells; endothelial cells [M3:20™]  |
| Skin (adult; newborn)  | Epidermal keratinocytes co-cultures; Dermal Fibroblasts; Mesenchymal cells [M3:10™]; others  |
| Tumors; various  | Rat, mouse, hamster; epithelial, mesenchymal, lymphoid; [M3:10™]; etc.   |

Legend to Tables 2 and 3. Supplemented M3 media formulations may contain various percentages of FBS added to M3Base™. Those media are designated as M3:{% serum}. As an example, M3:10 contains 10% v/v FBS. Some media are kept in stock, while others are special manufactured or made by the customers by adding the FBS to the media.

## Storage of M3:10AC™ Media

M3:10AC™ is refrigerated at 2°C to 8°C when not in use. Do not freeze. The shelf-life is 18 months from the manufacturing date.

## Master Files Applications Note

The M3™ Family of media is in FDA Drug and Device Master Files but have not been tested by INCELL for any specific diagnostic or therapeutic use. To request use of a Master File call, FAX, or email to [masterfiles@incell.com](mailto:masterfiles@incell.com).

## Ordering: Contact Charter Medical

Toll Free: 866.458.3116  
incell@chartermedical.com  
FAX: 336.714.4241

## Technical Assistance

The scientists at INCELL are available to discuss the media or special needs of your cells, and to assist you in your cell culture applications. Call 1-800-364-1765 or e-mail [info@incell.com](mailto:info@incell.com).